

- Unpack carefully the balance because it is a precision instrument.
- Put the balance on its place, which logically should be firm and horizontal.
- Place the pans so that the adhesive coincides with the one placed on the cross.
- To weigh with this type of balance, put the object to be weighed on the left pan. Immediately after compensate the right pan with weights.
- Only in the case that the balance loses its balance and in order to set it to zero, slide the counterpoise situated below.
- We grant a WARRANTY of 10 years for the balance model K-2000.
- The warranty cannot be awarded for damages caused by fall or bad use.

The balance is a precision instrument which has to be handled carefully in order to obtain the best performance during many years.

zu beziehen bei

sold by

www.conatex.com

