# QUICK START GUIDE

## EN

### Quick Start Guide

Thank you for purchasing this Moticam. This camera is a delicate instrument and should be handled with care. When not in use, please store the camera in a safe place with the dust cap attached. Do not let the camera get in contact with water or expose it to extreme temperatures (below 0 or above 50 degrees Celsius). In order to receive the highest level of service, please retain your receipt as proof of purchase and **register your camera online at [www.motic.com](http://www.motic.com) today.**

Should you have any questions please do not hesitate to **contact our customer service.**

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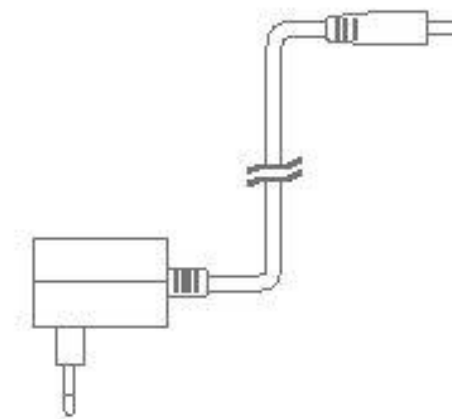
#### EN Using the camera on an eyepiece

1. Install the Motic Images Plus application program from the CD included.
2. Remove the Dust Cap (B) from the camera (A) and attach the focusable lens (C).
3. Choose the right size eyepiece adapter (D) and attach this to the focusable lens.
4. Attach this assembly to your microscope's eyepiece by tightening the centering screws on the eyepiece adapter.
5. Connect the USB cable (F) onto the camera and then plug it into the computer.
6. Install the driver if necessary by following the instructions on screen. The driver for the Moticam 580 is located in the Moticam 580 Subfolder on the CD.

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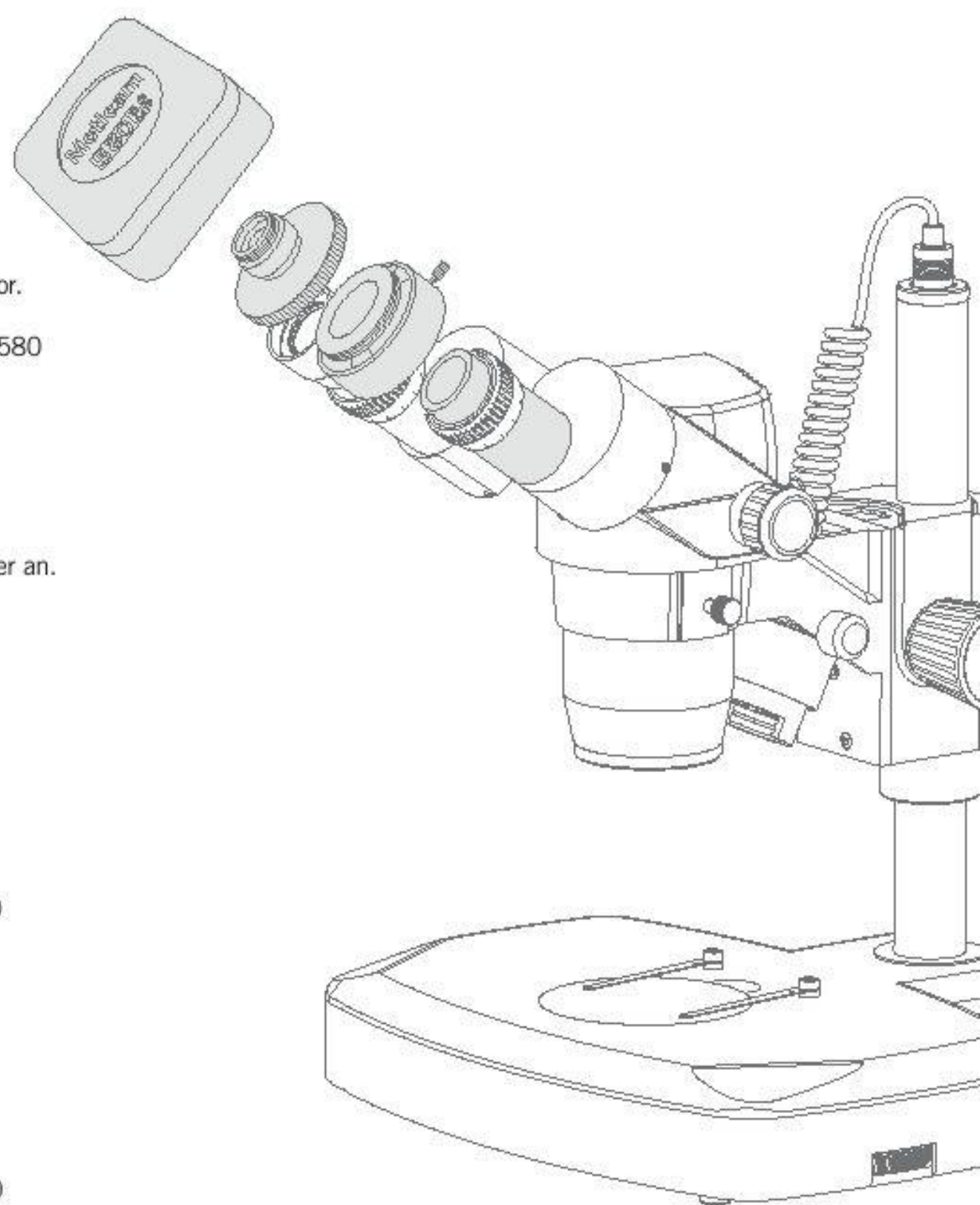
#### EN Using the camera as a Macro Camera

1. Install the Motic Images Plus application program from the CD included.
2. Remove the Dust Cap (B) from the camera (A) and attach the focusable lens (C).
3. Attach the Macro Tube (H) onto the lens.
4. Place this assembly over the item you wish to view and ensure there is adequate lighting. If necessary, a desk lamp is a good solution.
5. Connect the USB cable (F) onto the camera and then plug it into the computer.
6. Install the driver if necessary by following the instructions on screen. The driver for the Moticam 580 is located in the Moticam 580 Subfolder on the CD.

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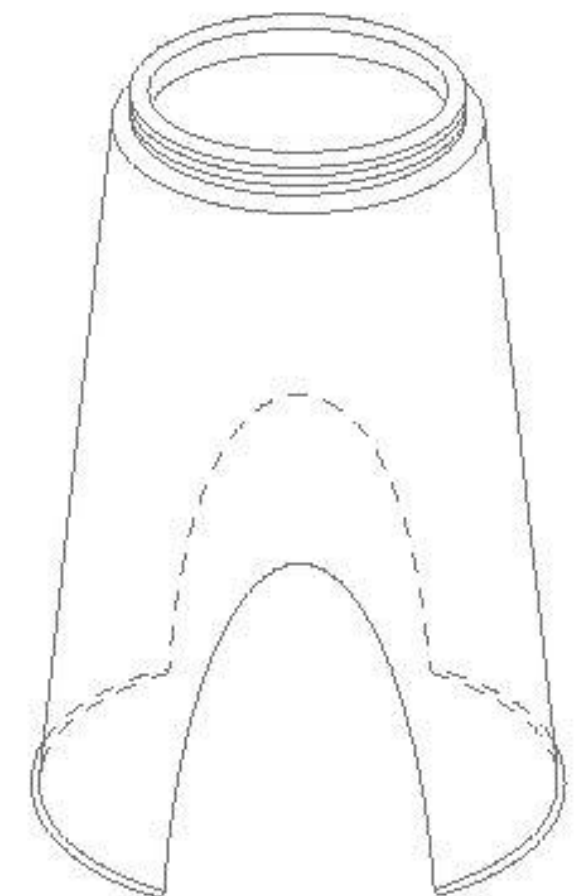
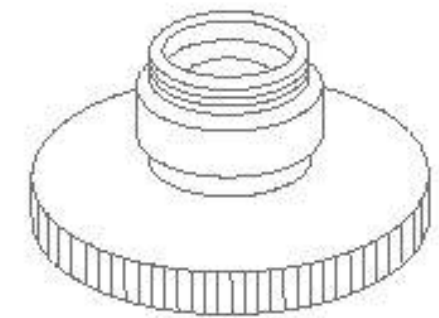
#### EN Using the camera on a Trinocular Microscope

1. Install the Motic Images Plus application program from the CD included.
2. Remove the Dust Cap (B) from the camera (A) and attach the microscope's C-Mount adapter (\*Please note that the C-Mount adapter is specific to your microscope. For Motic microscopes, please consult the relevant Motic Accessory list and choose an appropriate adapter for your Moticam. Depending on the type of adapter and microscope, the C-Ring (E) may be necessary between the camera and the C-mount adapter).
3. Connect the USB cable (F) onto the camera and then plug it into the computer.
4. Install the driver if necessary by following the instructions on screen. The driver for the Moticam 580 is located in the Moticam 580 Subfolder on the CD.

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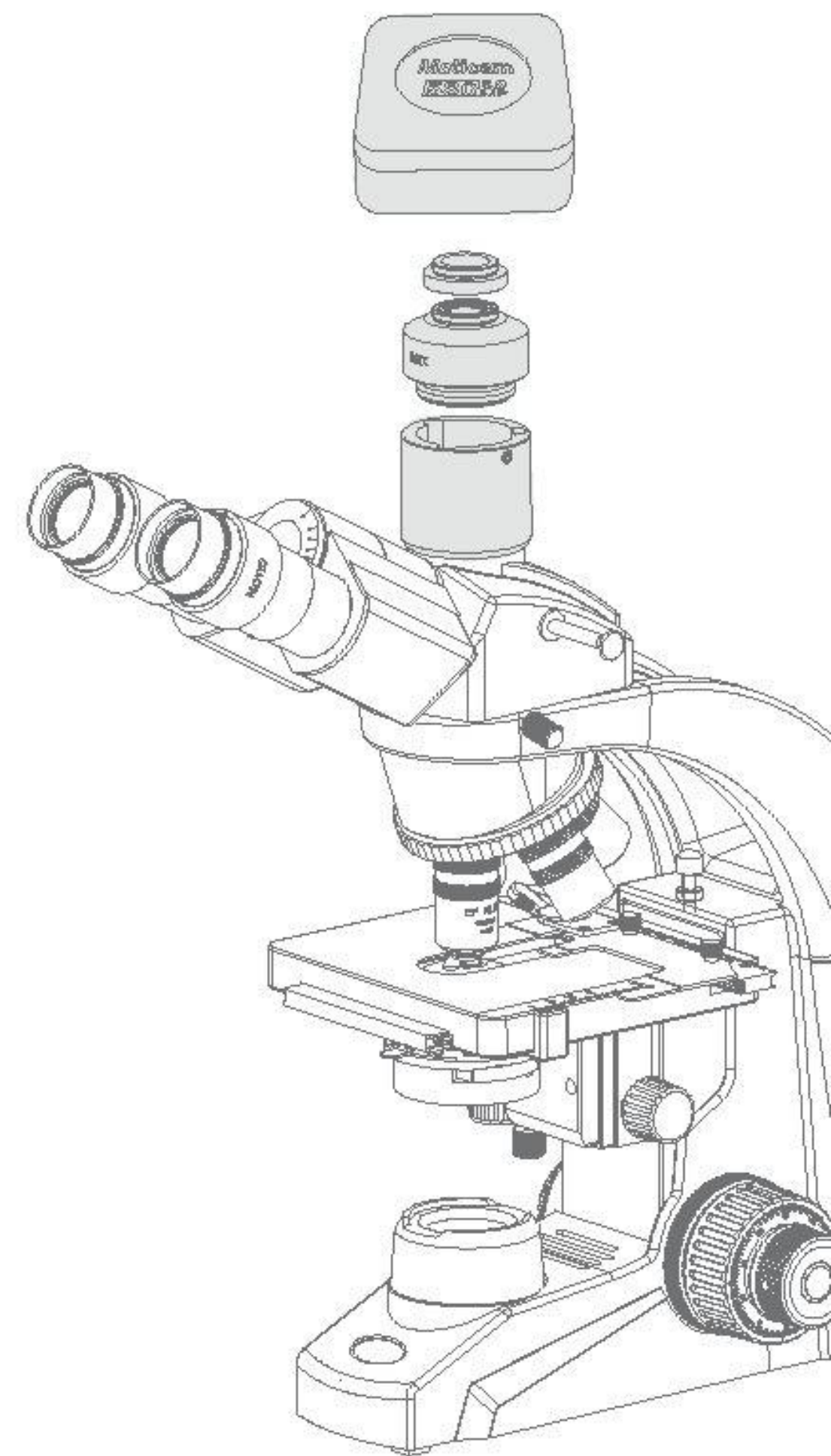
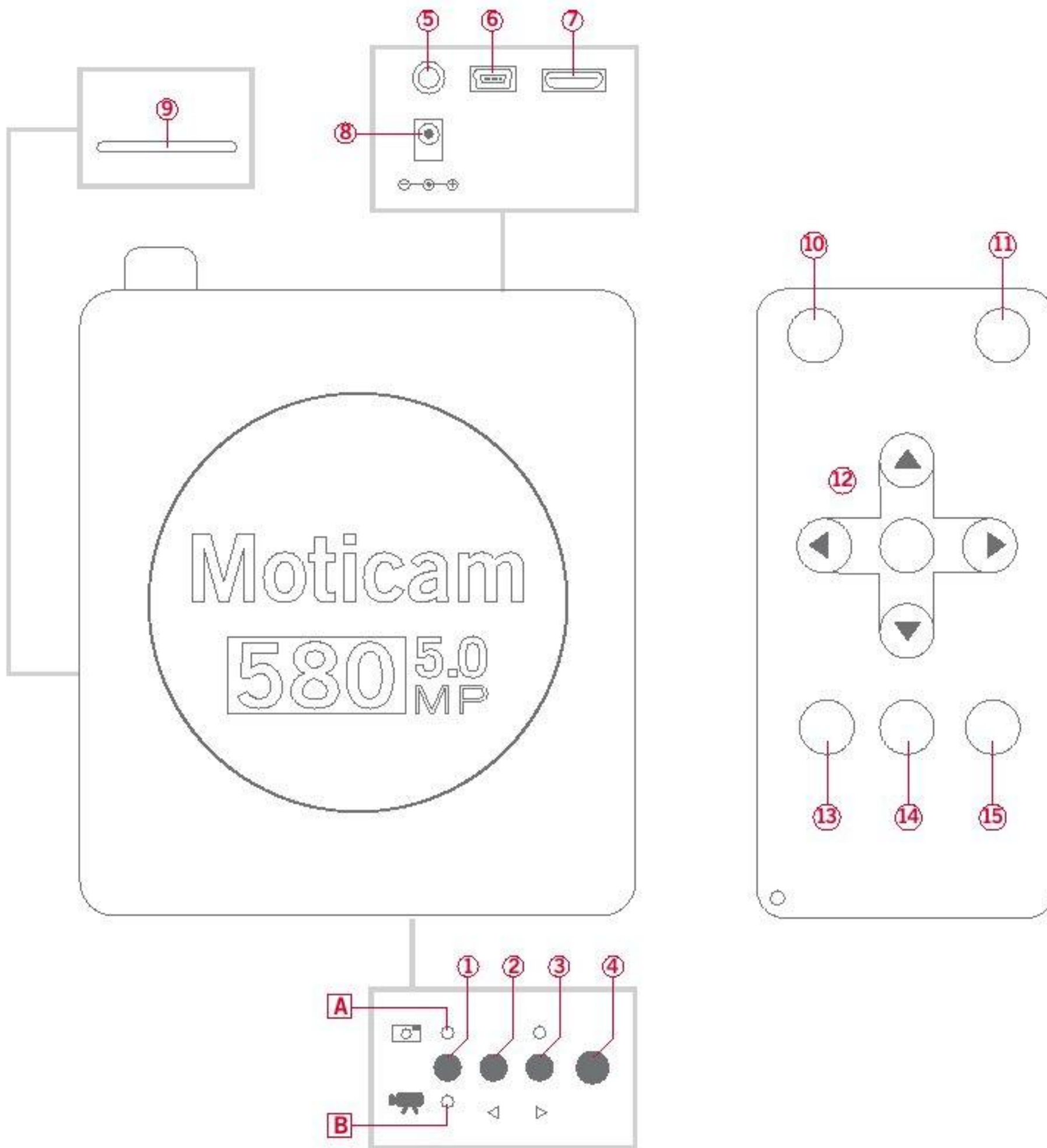
**EN** The button controls on the camera are only active when the Moticam 580 is connected to a Monitor using the HDMI or AV out connections. The Power Supply must be used in this case. When the camera is connected to a computer using the USB output, the Moticam 580 switches to PC Camera Mode and the camera can then be controlled through Motic Images Plus. When using the camera through a USB cable, no Power Supply is needed.

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**EN** 1. This button cycles through the functions of the camera. The LED indicator lights show what mode the camera is currently in. If Indicator A is on, the camera is in the Still Image Capture Mode. Pressing the SNAP button captures an image onto the SD Card. If Indicator B is on, the camera is in the Live Image Capture Mode. Pressing the SNAP button captures a video file onto the SD Card. If both Indicators A and B are on, the camera is in the ALBUM mode, use buttons 2 and 3 to scroll through the images on the SD Card.

2. When an analog device is connected to the camera, this button switches the output between NTSC and PAL outputs. When the camera is in ALBUM Mode use this button to scroll left.

3. Pressing this button activates the camera's hardware White Balance function. Use this function to calibrate a white background depending on the microscope's illumination. When the camera is in ALBUM Mode use this button to scroll right.

4. This button will Capture a Still Image or Start/Stop the Video Recording depending on what Mode the camera is in.

5. AV out for connecting to an Analog Display (the cable is provided in the Moticam 580 box)

6. USB port for connecting to a Computer (the cable is provided in the Moticam 580 box)

7. Mini-HDMI port for connecting to an HDMI ready Display (the cable is provided in the Moticam 580 box)

8. Power port for connecting the included Power Supply.

9. SD Card slot for capturing images and video files onto an SD card. The maximum size compatible is 32GB SDHC.

*NOTE: The remote control for the Moticam 580 is an optional extra available in the premium pack*

10. Pressing this button switches off the camera completely. Please note that you will need to replug the power supply to switch it on.

11. This button will Capture a Still Image or Start/Stop the Video Recording depending on what Mode the camera is in.

12. This toggle pad allows you to scroll through the images in the ALBUM and also through the camera's custom setup MENU.

13. Access the camera's custom setup by pressing this button. Please note that all settings go back to the default once the camera is disconnected from the power supply.

14. This button cycles through the functions of the camera. The LED indicator lights show what mode the camera is currently in. If Indicator A is on, the camera is in the

Still Image Capture Mode. Pressing the SNAP button captures an image onto the SD Card. If Indicator B is on, the camera is in the Live Image Capture Mode. Pressing the SNAP button captures a video file onto the SD Card. If both Indicators A and B are on, the camera is in the ALBUM mode, use buttons 2 and 3 to scroll through the images on the SD Card.

15. Pressing this button activates the camera's hardware White Balance function. Use this function to calibrate a white background depending on the microscope's illumination.

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